


How-to Instructions

BarbLock® Ultra-Secure Retainers

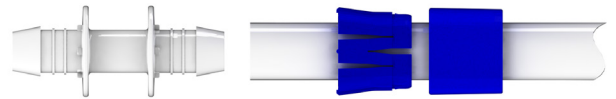


- 01 Load sleeve and collet on tube, leading with ends marked "Top".




Standard two-piece retainer


- 02 Insert barbed fitting into tube. Leave a space of approximately 1/16" between the tube and tube stop. This corresponds to the last fitting's rib.

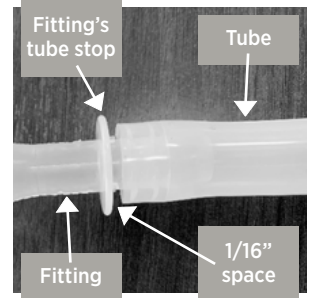


- 03 Move collet to base of tube. Leave the same approximately 1/16" space between collet and tube stop.

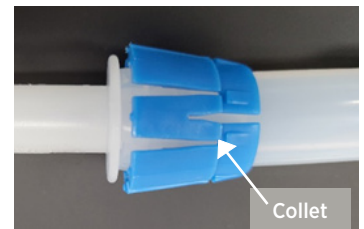

- 04 Lower sleeve over collet.

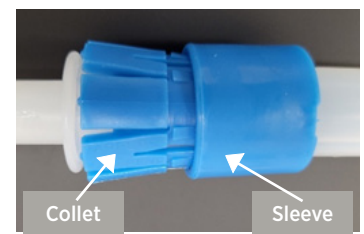

- 05 Lock assembly with a BarbLock® tool. The jaws should fit between the collet and tube stop, and behind the sleeve.

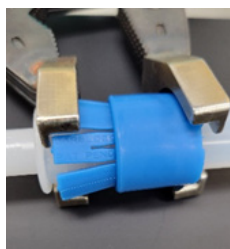

- 06 When the assembly is properly locked, tubing cannot be pulled. The space remains between the BarbLock® and tube stop, and tubing may expand in this space, but it should not be swollen.

